

Nr: 0012-R01e
2020-03-25

Operating Instructions

According to the
Betriebssicherheitsverordnung

UNIVERSITÄT
DUISBURG
ESSEN

Offen im Denken

Faculty of Chemistry

Institute: AAC

Work place: AAC Labs

Activity: Laboratory work

EQUIPMENT

Laboratory dishwasher Miele Professional g 7883

DANGERS FOR HUMANS AND ENVIRONMENT



- Danger due to glass breakage.
- Risk of burns due to alkaline detergent and citric acid.
- Residues of solvents can form explosive mixtures.
- Danger due to hot steam escaping when opening the machine after the rinsing process.

PROTECTIVE MEASURES AND RULES OF CONDUCT



- Wearing the PPE: gowns, glasses, protective gloves.
- Carefully sort glassware and watch out for any risk of injuries and contamination.
- Pre-clean glassware!
- Do not open the door during the flushing process!
- After the program has finished, the radiators inside are still hot for a few minutes, so do not touch them!
- Immediately after the end of the program, when the door is opened, hot steam will be released, so keep a safe distance.
- Do not remove glassware when hot, only in warm, better cold condition.

RESPONSE TO MALFUNCTIONS



Shut down device immediately, faults must be reported to:

- Maria Madani, if not present O. Schmitz, M. Sulkowski, F.Uteschil, S. Meckelmann.

BEHAVIOR IN CASE OF ACCIDENT / FIRST AID



- Keep calm.
- Call first responders.
- Emergency call: 0112
- Report accident.

INSTANDHALTUNG / ENTSORGUNG

- Maintenance only by authorized, competent persons.
- Switch off the device and disconnect it from the mains.
- Regular inspection of wearing parts.